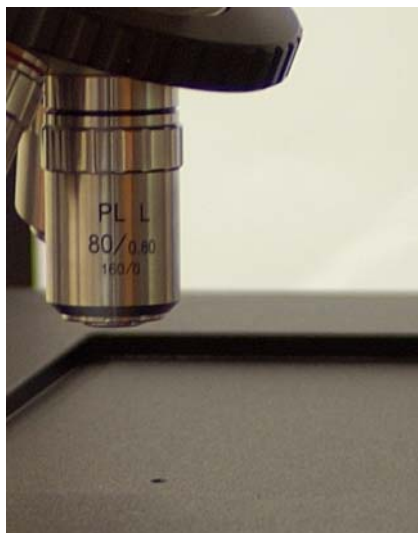
Microscopio Metalográfico, modelo 101 | Metallurgical microscope, model 101

Marca|Trademark:Zuzi

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Microscopio Metalográfico, modelo 101 | Metallurgical microscope, model 101

Marca|Trademark:Zuzi

Zuzi metallurgical microscope model 101 has been especially conceived for observation of any kind of opaque surfaces such as metals, alloys or mineral, by means of the incidence of a light beam that is reflected by the sample and reaches the objectives. Thanks to its excellent technical features it is a equipment with a wide range of applications in industry, materials science, geology, etc.

Ergonomics

It is a very solid, large and stable equipment as well as with a comfortable and useful design. The head is 30° inclined and presents interpupillary distance adjustment that together with the diopter correction ring in one of the eyepieces allows an adequate observation. It is provided with a third tube for graphically recording images of the samples by using a digital photograph or video camera thus increasing considerably the equipment features. The light beam is directed by means of a sliding rod that sends 100% of the light to the eyepieces or to the third tube.

Sample focusing is performed through coarse and fine focusing knobs placed at both sides of the microscope in a lower position for a more comfortable handling. Moreover, the right knob is provided with a friction control ring and the left one with a focusing upper limit blocking lever to avoid the slide to accidentally hit the objectives.

Triple layer stage

The mechanical stage is very large (250x230 mm) and can be moved 154 mm along both axes. A special mechanism is used for moving the sample: the triple layer stage. It presents two movements in common with the double layer stages, but it is also provided with a third layer specialized in quick movements all along the stage, very useful for studying large samples.

Illumination

Zuzi metallurgical microscope model 101 is equipped with a Köhler illumination system that allows a better control over the light that reaches the sample. This illumination system includes a halogen lamp with intensity adjustment, a horizontally adjustable condenser and field and aperture diaphragms. Moreover, the lamp is provided with a centering mechanism to get a correct adjustment and thus assuring an optimum illumination and visualization of samples. The microscope also presents a filter holder with green, yellow, blue and matt glass filters to obtain a higher quality image.

Polarization

The microscope also allows the study of the optical properties of samples since it is equipped with an analyzer and a polarizer that can be rotated until polarizer and analyzer are orthogonally positioned. Both, analyzer and polarizer can be removed from the light path when they are not being used.

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